



RCPN

Standard operating procedure of Micro centrifuge

Objective: This instruction explains how use and keep the micro centrifuge apparatus.

Scope: using for centrifuging the different ranges of solutions with maximum 2 mL volume.

Responsibility: Laboratory assistance is responsible for the accuracy of the apparatus and should be reported to him/her in case of problems.

Materials and equipment: ---

Guidelines for use:

- 1- Turn on the apparatus.
- 2- Open the door of apparatus with blue key.
- 3- Place the tubes in face to face pattern for maintaining the balance of apparatus.
- 4- Adjustment the time and centrifuging speed of the apparatus with the blue set bottoms and run your reaction.

Warning and safety precautions:

If any abnormal sound or vibrate are felt, press the “Stop” bottom immediately.

Maintenance:

The door of the apparatus should be always closed.

Document ID: **RCPN 43**

Document name	Standard operating procedure of Micro centrifuge
Document arrangement	Mohanna Osali
Date	10/25/2018
Full file name	Standard operating procedure of Micro centrifuge
Document Description	This instruction explains how use and keep the micro centrifuge apparatus.
Author / translator	Rahimeh mousavi